



BioClean-D™ Sterile Disposable Coverall with Collar DONNING PROCEDURE



1 Don first pair of BioClean™ sterile gloves, facemask and bouffant cap. Don sterile BioClean-D hood.

2 Open vacuum-packed bag using EasyTear™ edge. Remove inner pack and open using EasyTear™ edge.

3 Remove coverall from inner pack, gently unfold and unzip ensuring it does not touch the floor, before holding the coverall inside at the waist.



4 Put one foot through ankle opening, pull coverall up over leg. Do the same with the other foot.

5 Insert one hand through wrist opening, bringing coverall up over arm. Do the same with the other arm. Put thumbs through the thumb loops.

6 Cross legs, straighten back and zip up coverall.

STEP
— by —
STEP



7 Remove backing tape on zip flap and press down lightly. Sanitize gloves.



8 Don sterile BioClean-D™ overboots using aseptic technique. Please refer to BioClean-D™ Sterile Overboots donning procedure.



9 If required don BioClean sleeve covers using aseptic technique. Please refer to BioClean-D™ Sterile Sleeve Covers donning procedure.



10 Don BioClean goggles.



11 Sanitize gloves. Don second pair of sterile gloves, using aseptic donning technique.



12 Perform a final inspection prior to entry into the sterile cleanroom, ensuring no skin is visible and all apparel is donned correctly.