



# BioClean-D™ Sterile Disposable Coverall with Hood DONNING PROCEDURE



**1** Don first pair of BioClean™ sterile gloves, facemask and bouffant cap.



**2** Open vacuum-packed bag using EasyTear™ edge. Remove inner pack and open using EasyTear™ edge.



**3** Remove coverall from inner pack, gently unfold and unzip ensuring it does not touch the floor, before holding the coverall inside at the waist.



**4** Put one foot through ankle opening, pull coverall up over leg. Do the same with the other foot.



**5** Insert one hand through wrist opening, bringing coverall up over arm. Do the same with the other arm. Put thumbs through the thumb loops.



**6** Bring hood over head touching only the inside.

STEP  
—by—  
STEP



**7** Cross legs, straighten back and zip up coverall.



**8** Remove backing tape on zip flap and press down lightly. Sanitize gloves.



**9** Don sterile BioClean-D™ overboots using aseptic technique. Please refer to BioClean-D™ Sterile Overboots donning procedure.



**10** If required don BioClean sleeve covers using aseptic technique. Please refer to BioClean-D™ Sterile Sleeve Covers donning procedure.



**11** Don BioClean goggles.



**12** Sanitize gloves. Don second pair of sterile gloves, using aseptic donning technique.



**13** Perform a final inspection prior to entry into the sterile cleanroom, ensuring no skin is visible and all apparel is donned correctly.