



BioClean-D™ Sterile Drop-Down Garment DONNING PROCEDURE




1 Remove drop-down coverall.



2 Hold internal red tab in your right hand and white tab in your left. Shake the garment to un-fold.



3 Insert one arm and then the other. Put thumbs through thumbloops

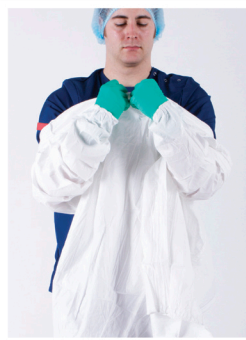


4 Hold the inside of the coverall and bring over your head.



Shake garment down allowing it to drop-down over body or use external tabs to pull garment down.



5 Put right foot through ankle opening and then foot-loop.

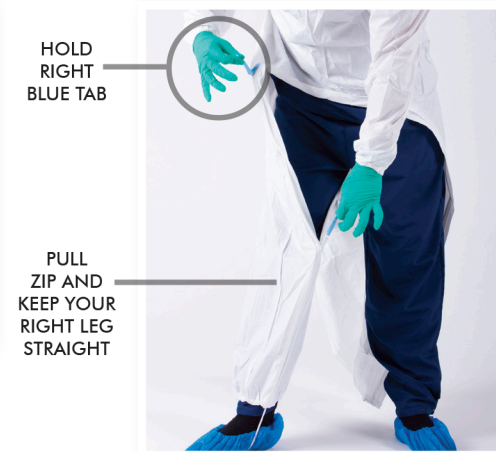

6 Hold blue tab on right side of waist. Pull zip up ensuring you keep your right leg straight.



7 Still holding blue tab pull zip round to the blue tab on the left hand side of waist.



8 Pull off blue tab on the right hand side.



9 Hold the blue tab on the left side of waist with your left hand.



Pull zip down with your right hand and remove the blue tab at the waist as you do so.



10 Remove the zip tab by pulling the tab through the zip hole. Discard all tabs.



11

Don sterile BioClean-D[™] overboots using aseptic technique. Complete gowning by donning goggles and a second pair of sterile gloves.